

UV-VIS LIGHT SOURCE

HIGH POWER UV-VIS FIBER LIGHT SOURCE

L10290

OVERVIEW

The L10290 is a newly developed UV-visible fiber light source that outputs 200 nm to 1600 nm light through a light guide (sold separately).

Using a "high brightness lamp", the L10290 provides twice radiant intensity that of conventional models. Besides a compact size and light weight that make it simple to carry, the L10290 is specifically designed for easy use. These features make the L10290 ideal for all types of portable devices.



* Light guide is sold separately.


FEATURES

- High Output: Twice Intensity (Compared to Conventional Model)
- High Stability: Fluctuation 0.004 %p-p Typ. (Equivalent to 2×10^{-5} A.U.)
- Long Life Lamp: 2000 Hours
- External Control
- Shutter Function
- Filter Holder (Sold Separately)


APPLICATIONS

- Spectrophotometry
- Environmental Measurement
- Absorption Spectrum Measurement
- Microtiter Plate Reader
- Pharmaceutical Testing
- Film Thickness Measurement
- Semiconductor Inspection
- Optical Component Inspection

SPECTRAL DISTRIBUTION (TYPICAL DATA)

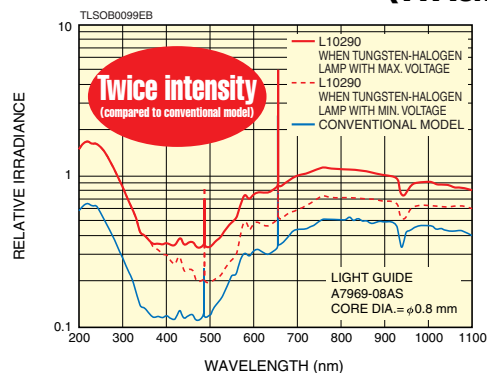

LIGHT OUTPUT STABILITY (at 230 nm)

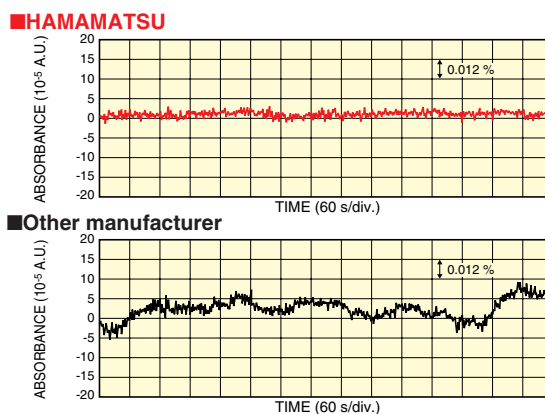

TLSOB0059EB

SPECIFICATIONS

Parameter		Specifications
Spectral Distribution		200 nm to 1600 nm ①
Built-in Lamp		High Brightness Deuterium Lamp 9 W Tungsten Halogen Lamp ②
Output	Fluctuation (Typ.)	0.004 % p-p (Equivalent to 2×10^{-5} A.U.)
Stability	Drift (Max.)	± 0.3 %/h
Input Voltage (AC)		100 V to 240 V (100 V/200 V Auto Switching) Single Phase 50 Hz/60 Hz
Power Consumption (Max.)		140 VA
Weight		Approx. 4 kg
Light Guide Connector		SMA905D
Shutter		Included
External Control		Lamp ON/OFF ③
		Shutter open/close
		Status Signal for Lamp ON/OFF and Main Power ON/OFF
Main Unit Switch		Main Power ON/OFF
		Lamp ON/OFF ③
		Shutter open/close
Applicable Standard		CE Marking


NOTE: ① The deuterium lamp (200 nm to 400 nm) and tungsten halogen lamp (400 nm to 1600 nm) can each radiate light separately.
 ② Tungsten halogen lamp intensity is adjustable within a range of ± 20 %.
 ③ The deuterium lamp and tungsten halogen lamp can each be turned on or off separately.

REPLACEMENT LAMP

Type No.	Lamp	Lamp Life ④
L10295	High Brightness Deuterium Lamp	2000 h
L10296	9 W Tungsten Halogen Lamp	2000 h

NOTE: ④ Guaranteed life. Defined as the time when the radiant intensity falls to 50 % of its initial value.

DIMENSIONAL OUTLINES (Unit: mm)

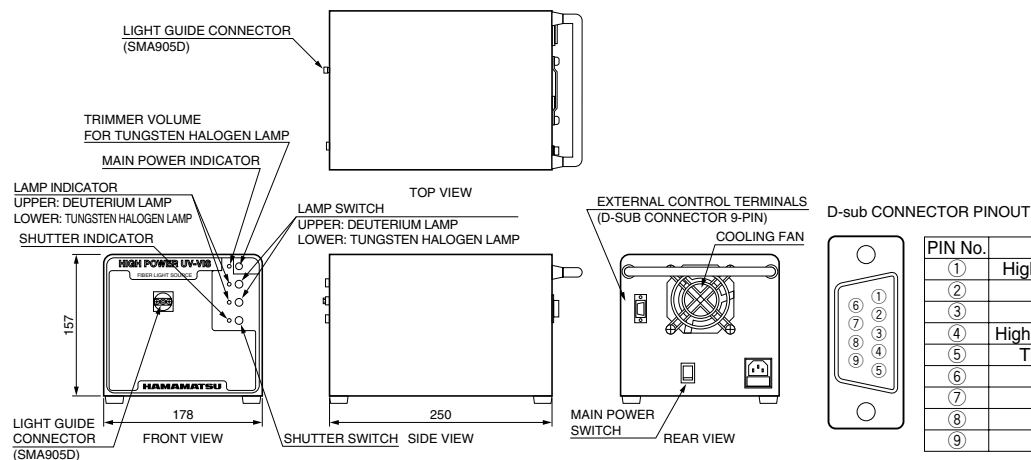

CE This product complies with safety standard EN61010-1, EMC standard EN61326-01.

Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult with our sales office.
 Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications are subject to change without notice. No patent rights are granted to any of the circuits described herein. ©2011 Hamamatsu Photonics K.K.

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P. O. Box 6910, Bridgewater, N.J. 08807-0910, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-2658 E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trappu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: infos@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road Welwyn Garden City Hertfordshire AL7 1BW, United Kingdom, Telephone: 44-(0)1707-294888, Fax: 44-(0)1707-325777 E-mail: info@hamamatsu.co.uk

North Europe: Hamamatsu Photonics Norden AB: Smidesvägen 12, SE-171-41 SOLNA, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01 E-mail: info@hamamatsu.se

Italy: Hamamatsu Photonics Italia: S.R.L.: Strada della Moia, 1/E, 20020 Arese, (Milano), Italy, Telephone: (39)02-935 81 733, Fax: (39)02-935 81 741 E-mail: info@hamamatsu.it


LIGHT GUIDE^⑤ (SOLD SEPARATELY)

Type No.	Core Diameter	Length
A7969-02AS ⑥	0.2 mm	1.0 m
A7969-06AS ⑥	0.6 mm	
A7969-08AS ⑥	0.8 mm	

NOTE: ⑤ Other shapes and combinations are available upon request.

⑥ UV resistant light guide.

UV RESISTANT CHARACTERISTICS (at 215 nm)

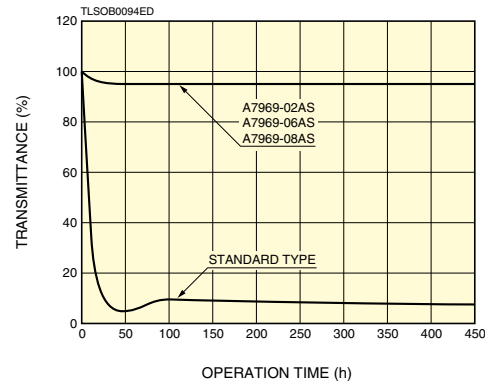

FILTER HOLDER (SOLD SEPARATELY)

Type No.	Applicable Filter
A10394	3 mm thick, 1 inch dia. or 1 × 1 inch

Filters should be prepared by the user.

When attaching a filter holder, remove the cover and insert the filter holder into the lamp house.

PIN No.	Signal Name
①	High brightness deuterium lamp ON/OFF
②	Tungsten halogen lamp ON/OFF
③	Shutter open/close
④	High brightness deuterium lamp status signal
⑤	Tungsten halogen lamp status signal
⑥	Main power ON/OFF status signal
⑦	No connection
⑧	No connection
⑨	GND

TLSOA0092EB

TLSZ1010E01
JUL. 2011 IP